FT 2

Single or dual drum tester with calculation of percent friability / weight loss. Print test protocols and connect your external analytical balance.



Friability and Abrasion Tester

FT2 — Single or dual drum testing.

sotax.com/FT2

Calculation of percent friability / weight loss and printing of test protocols

Perform two friability / abrasion tests simultaneously

Programmable rotational speed / time

Menu-guided qualification procedure with printed report

Integrated 10° tilting mechanism as per USP

Collection tray for each drum



User-friendly SingleButton™ navigation

Networking and data management with q-doc®

1-2 drums

Connect analytical balance

Print test protocols

Password protection

Electronic data management

Recording of weight readings from external analytical balance



Ease of regulatory compliance.



Prevent errors.


Manual tasks typically associated with the recording of weight measurements or report creation can be virtually eliminated by connecting a standard laboratory balance. If desired, operators can start tests using a predefined USP/Ph.Eur. method – ensuring that tests are performed according to specifications.

Robust design.

The FT2 features a maintenance-free drive that guarantees gentle starting and stopping motions. A quick-lock system prevents any slippage. Whether operated in single or dual configuration, friability drums don't have to be removed to insert tablets. On completion of a test run, samples are automatically discharged into separate collection trays.

Protocol your test results.

Test protocols including weight loss and percent friability calculations can be conveniently printed directly from the tester. Additionally, the qualification procedure is fully menu-guided including a printed report. If needed, the FT2 can also be integrated with q-doc® data management for consolidated batch reporting of various tests performed on different instruments.

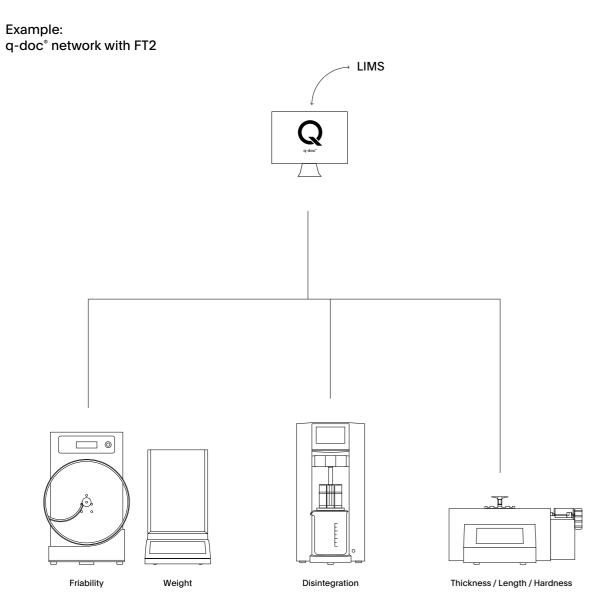

FT2 Technical Specifications

| Number of drums | | 1-2 |
|----------------------|---|----------------------------|
| Drum types | Friability drum, anti-static | 0 |
| | Abrasion drum, anti-static | 0 |
| Test modes | USP <1216> | • |
| | Ph. Eur. 2.9.7 | • |
| | User-defined | • |
| Rotation | Speed | 20 - 100 rpm, programmable |
| | Accuracy | ±1 rpm |
| Test modes | Time | • |
| | Revolution count | • |
| 10° tilting function | | • |
| Balance connection | Mettler Toledo* / Acculab* | • |
| User interface | Password protection | • |
| Printer | LAN printer / Epson* 40col. | Ethernet / RS-232 parallel |
| Data Management | q-doc*software, for data collection | 0 |
| Interfaces | | Ethernet, 2×RS-232 serial |
| Power supply | | 100 – 240 V / 50 – 60 Hz |
| Weight | (without packaging) | 8 kg (17.7 lbs) |
| Dimensions | Width | 305 mm (12 inch) |
| | Depth | 280 mm (11 inch) |
| | Height | 475 mm (18.7 inch) |
| CE Conformity | The FT2 fully complies with all CE and EMC equipment guidelines relative to mechanical and electrical safety and electromagnetic compatibility. | |

optional

quired Technical specifications are subject to change without prior notice. Products illustrated in this brochure may include options or modifications not fitted as standard. No liability for errors and omissions.

SOTAX Worldwide.

Europe

Switzerland (HQ Europe)

Aesch / Basel P Tech Support +41 61 487 5460 P Office +41 61 487 5454 info@sotax.com

Czech Republic

Prague P Tech Support +41 61 487 5460 P Office +420 246 039 260 sotaxCZ@sotax.com

France

Saint-Louis P Tech Support +41 61 487 5460 P Office +33 3 8970 0846 info@sotax.com

Germany

Lörrach P Tech Support +41 61 487 5460 P Office +49 7621 16 5635 info@sotax.com

Great Britain

Foston, Derbyshire P Tech Support +44 1332 981 571 P Office +44 1332 981 303 sotaxGB@sotax.com

Italy

Milan P Tech Support +41 61 487 5460 P Office +39 02 8363 2641 sotaxitalia@sotax.com

Ireland

Longford, Co. Longford P Tech Support +353 43 334 7779 P Office +353 43 334 7779 sotaxIE@sotax.com

Americas

USA (HQ Americas)

Westborough, MA
P Tech Support +1 508 544 4040
P Office +1 508 417 1112
sotaxusa@sotax.com

Canada

Westborough, MA
P Tech Support +1 508 544 4040
P Office +1 508 417 1112
sotaxcanada@sotax.com

Asia-Pacific

Thailand (HQ Asia-Pacific)

Bangkok P Tech Support +91 99 2021 1211 P Office +66 20 02 71 15 sotaxasia@sotax.com

India

Mumbai P Tech Support +91 99 2021 1211 P Office +91 22 2685 1903 sotaxindia@sotax.com